

-STR Thin Semrock filter installation instructions

1. You will need both filter wrenches pictured here. The standard filter wrench is visible on the right and the new filter wrench for the –STR filters on the left. **Remove the cover from the loading port**



2. Use the cup side of the standard filter wrench to **remove the cup from the filter wheel.**




3. With the **entire cup** removed from the filter wheel you should visibly see the filter wheel itself through the loading port.



4. Orient the new filter wrench vertically and **install the threaded filter on the wrench**. The design of the filter wrench should allow some flexibility to secure the threaded filter in place.





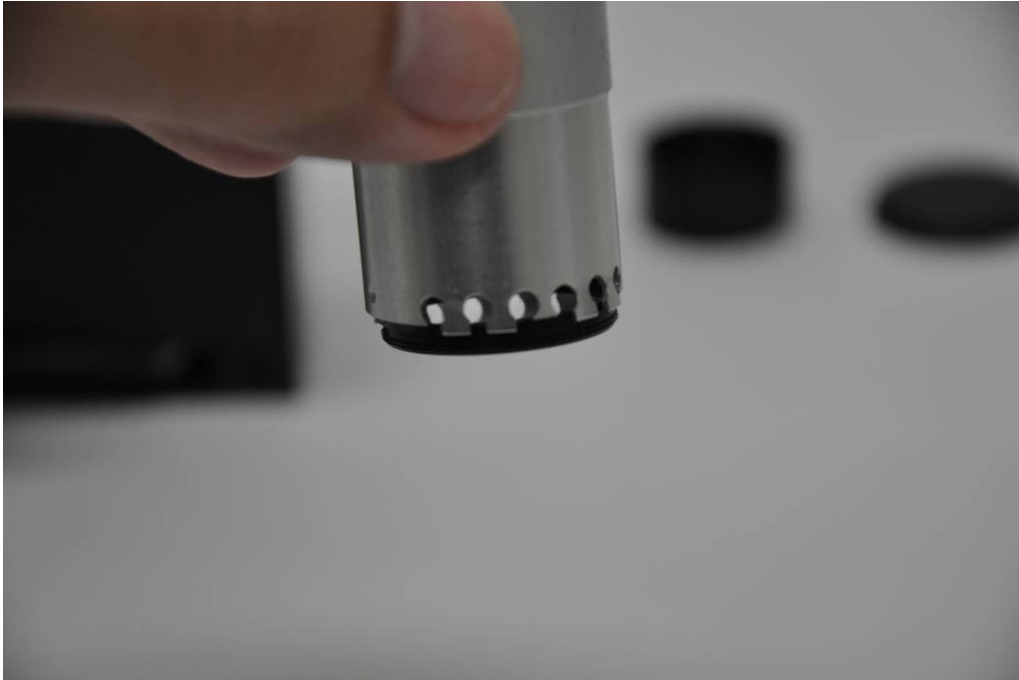

5. With the threaded filter secured to the wrench, **thread the filter directly into the filter wheel.**

